



Catalogue Number: AS-25241

Ninhydrin Test Kit for Resin

- **Description:** This kit consists of three bottles (10 ml each) of reagents. It is designed to check if the coupling of a protected amino acid to the amino group of the resin is complete.
- **Storage:** 2-8 °C
- **Test Protocol:**
 1. Place 20-30 beads of resin in a 12x75 test tube
 2. Add 2 drops of reagent 1, 2 and 3 to the test tube
 3. Heat the test tube in a heating block at approximately 120 °C for 3 minutes
 4. **Data Analysis**
 - 4.1 Colorless: Complete coupling
 - 4.2 Blue color: Incomplete coupling

AnaSpec Inc.

34801 Campus Drive, Fremont, CA 94555

Toll-Free: 800-452-5530 • Tel: 510-791-9560 • Fax: 510-791-9572

service@anaspec.com • www.anaspec.com